



ECD EFC INSTALLATION

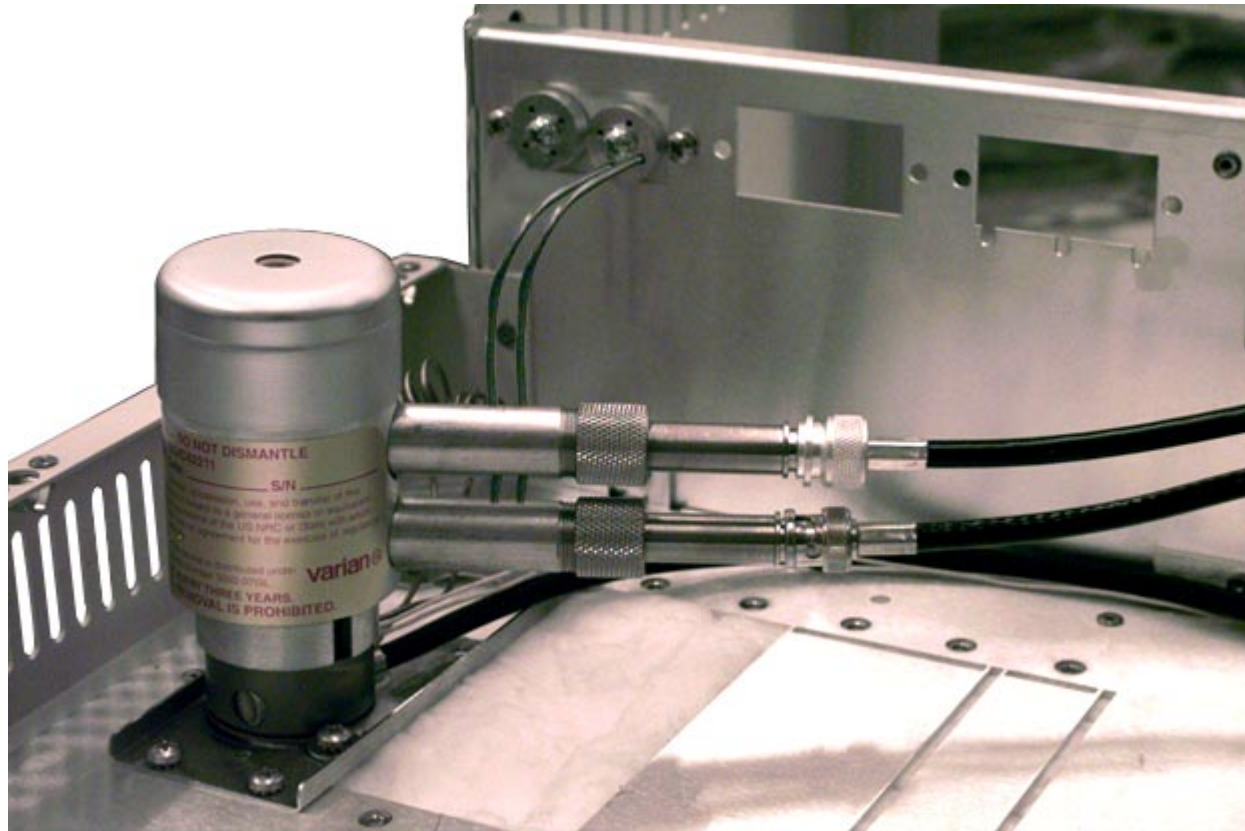

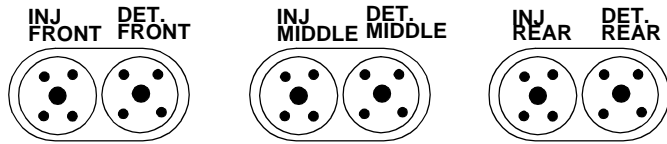


MANIFOLD PLUMBING

03-917142-00
VITON FERRULE


12-901150-00
SCREW

03-917141-00
CONDYNE PLATE


REF
ONLY

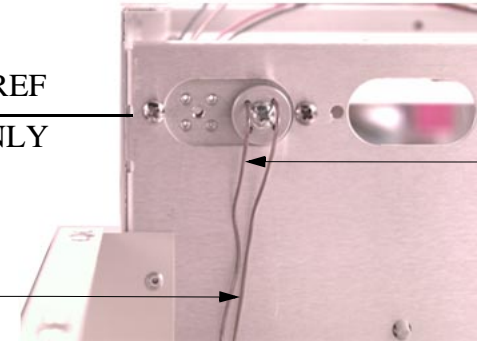

BOTTOM
LINE FROM
DET. BASE,
BLUE 

TOP
LINE FROM
DET. BASE,
(BLANK) 

DETECTOR SIDE MANIFOLD

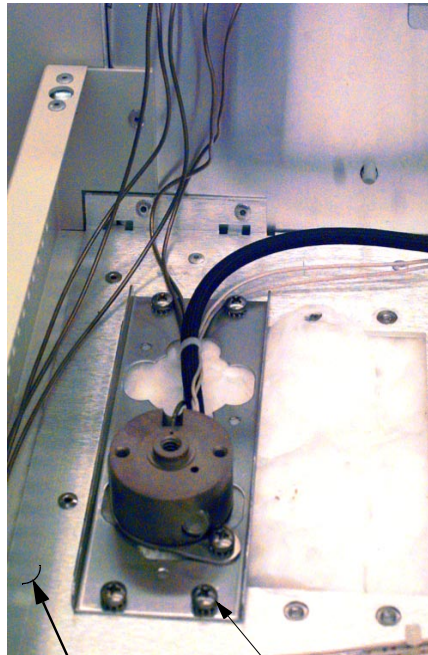

REF. OUTER
OVEN TOP


12-901163-00
SCREW, 4EA

03-925374-01
TEMPERATURE PROBE

03-925396-01
120V HTR. CART.

03-925396-02
230V HTR. CART.

03-925396-03
101V HTR. CART.

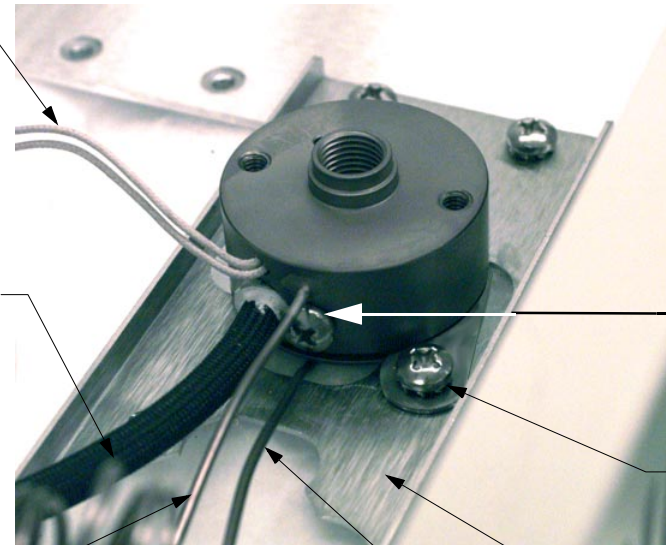

03-918663-06
SCREW

12-901163-00
SCREW, 2EA

03-925526-01
DETECTOR BASE CLIP





03-925326-01
DETECTOR MOUNTING BASE


DETECTOR BASE MOUNTING

[CLICK HERE FOR PROBE CONNECTION TO J7 ON MAIN BD.](#)

[CLICK HERE FOR HEATER CONNECTION J41 ON POWER BOARD](#)

CABLE INSTALLATION

CABLE PLUG-IN
TO MAIN BOARD

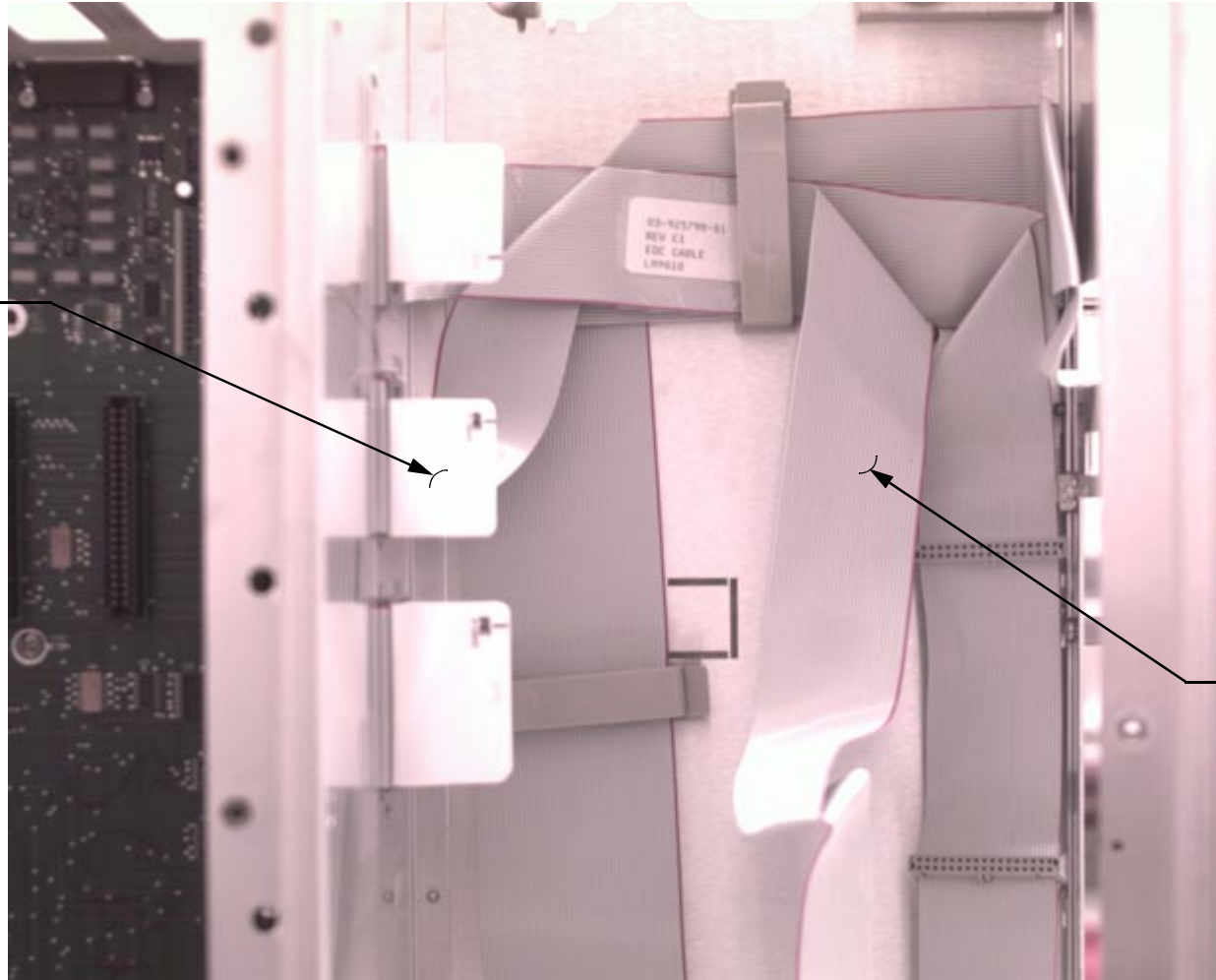

03-925798-01
CABLE, FLAT FOR
DETECTOR EFC


INLET MANIFOLD SUBASSEMBLY

03-925842-24
TUBING, PEEK 1/8 26" (NATURAL)


03-925838-23
TUBING, PEEK 23" (BLUE)

03-917084-00, 3 PLCS
NUT, VITON FERRULE
03-917157-00, 3 PLCS
WASHER, FERRULE
03-917142-00, 3 PLCS
VITON, FERRULE


03-918153-00
INLET MANIFOLD

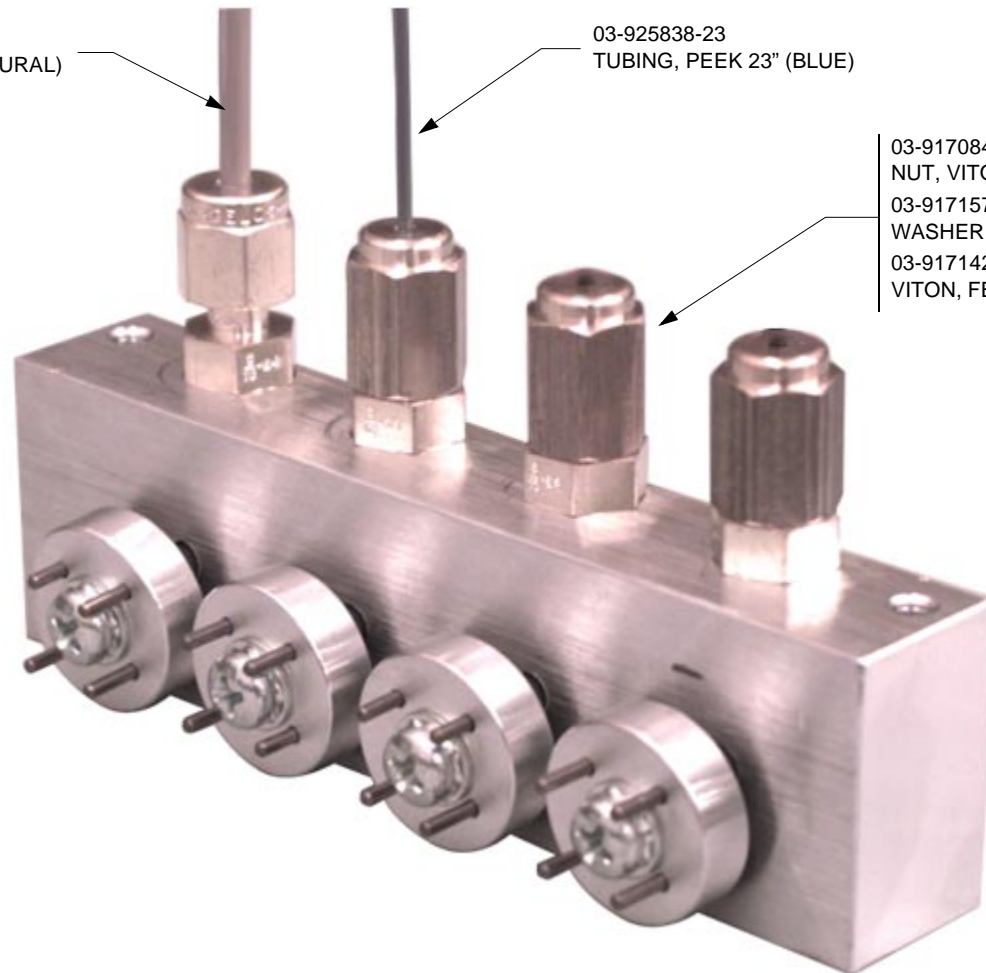

MODULE INSTALLATION

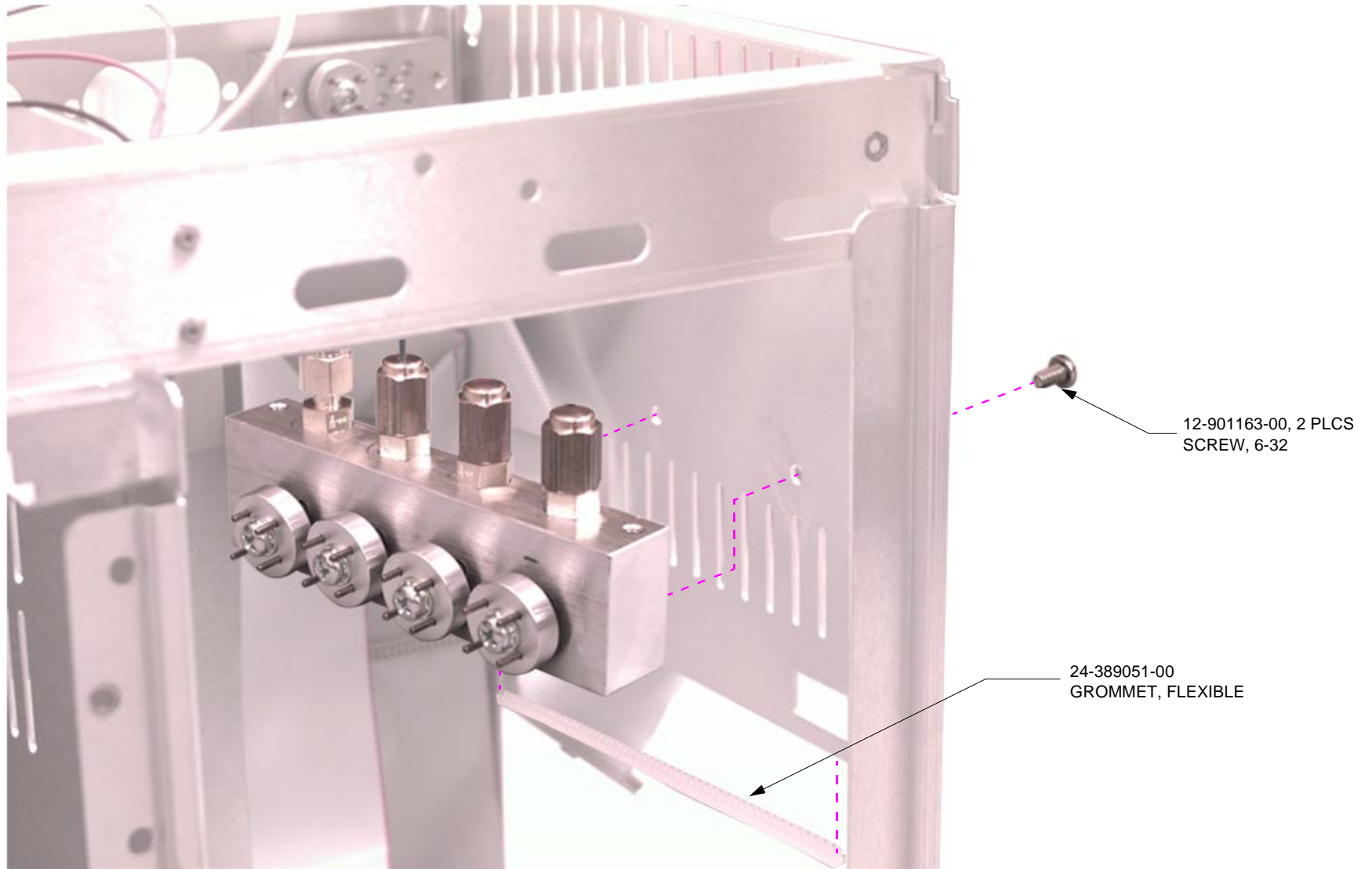


INLET MANIFOLD INSTALLATION

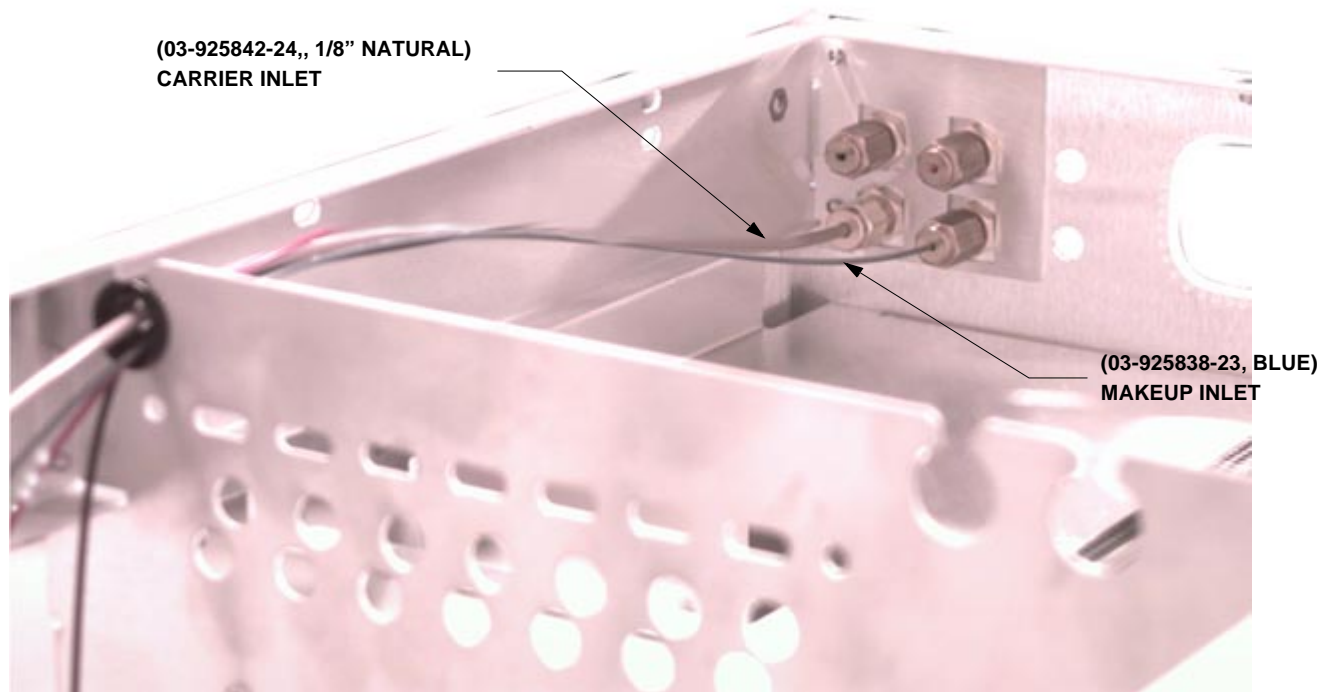

MODULE INSTALLATION

NOTES:

1. CONNECT CABLE TO MODULE, THEN MOUNT MODULE TO FRONT PANEL.
2. MOVE MODULE LOCATION JUMPER TO SPECIFY FRONT, MIDDLE, OR REAR.

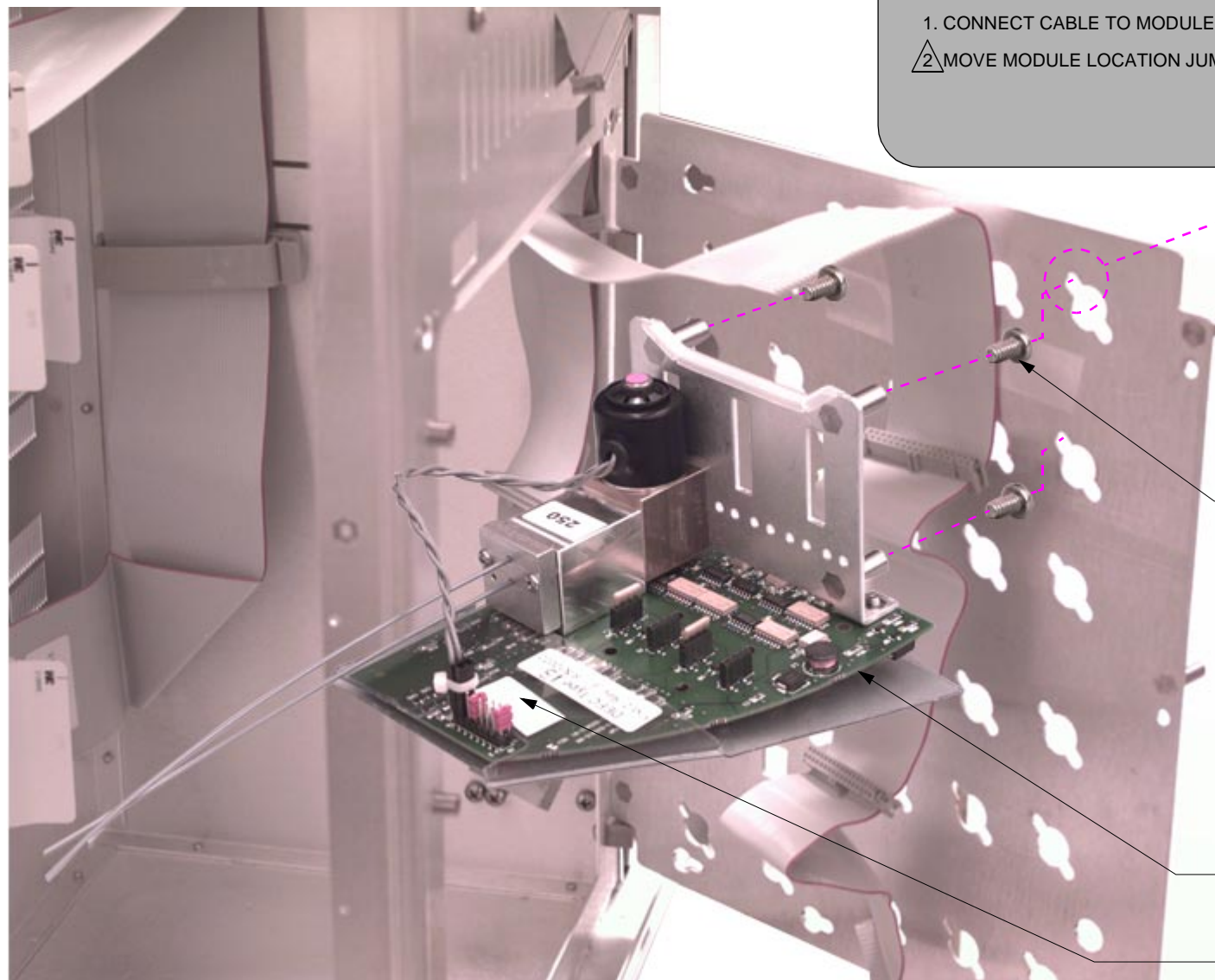


NOTE MOUNTING LOCATION

12-901163-00, 4 PLCS SCREW, 6-32

03-925609-01 EFC MODULE, FID


2

ADDITIONAL MODULE CABLEING (IF REQ'D)

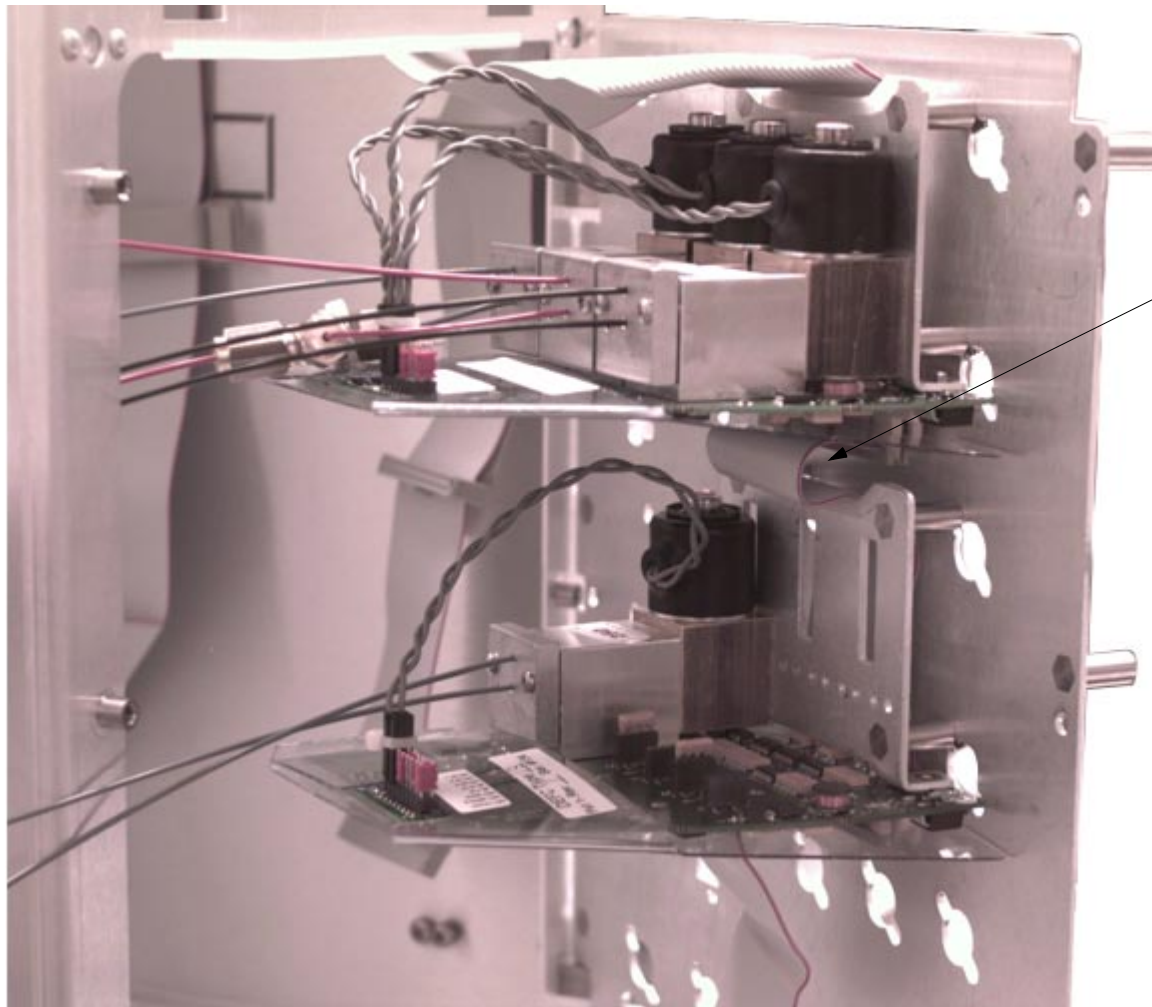

NOTE CABLE SERVICE
LOOP FOR ADDITIONAL
DETECTOR EFC MODULE

PNEUMATICS SIDE MANIFOLD

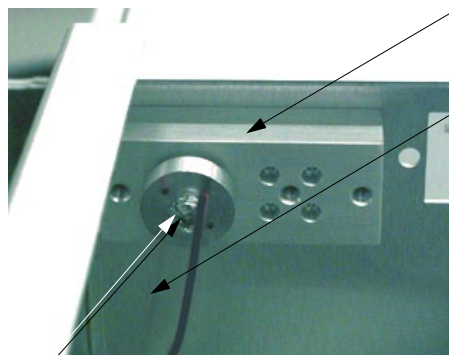

REFERENCE ONLY
PNEU. MANIFOLD

MAKE UP LINE
FROM MODULE TO
DETECTOR BLUE (26")

03-917142-00
VITON FERRULE


12-901150-00
SCREW

03-917141-00
CONDYNE PLATE

REF. ONLY


MODULE PNEUMATICS CONNECTIONS

MAKEUP LINE TO MODULE, BLUE (16")

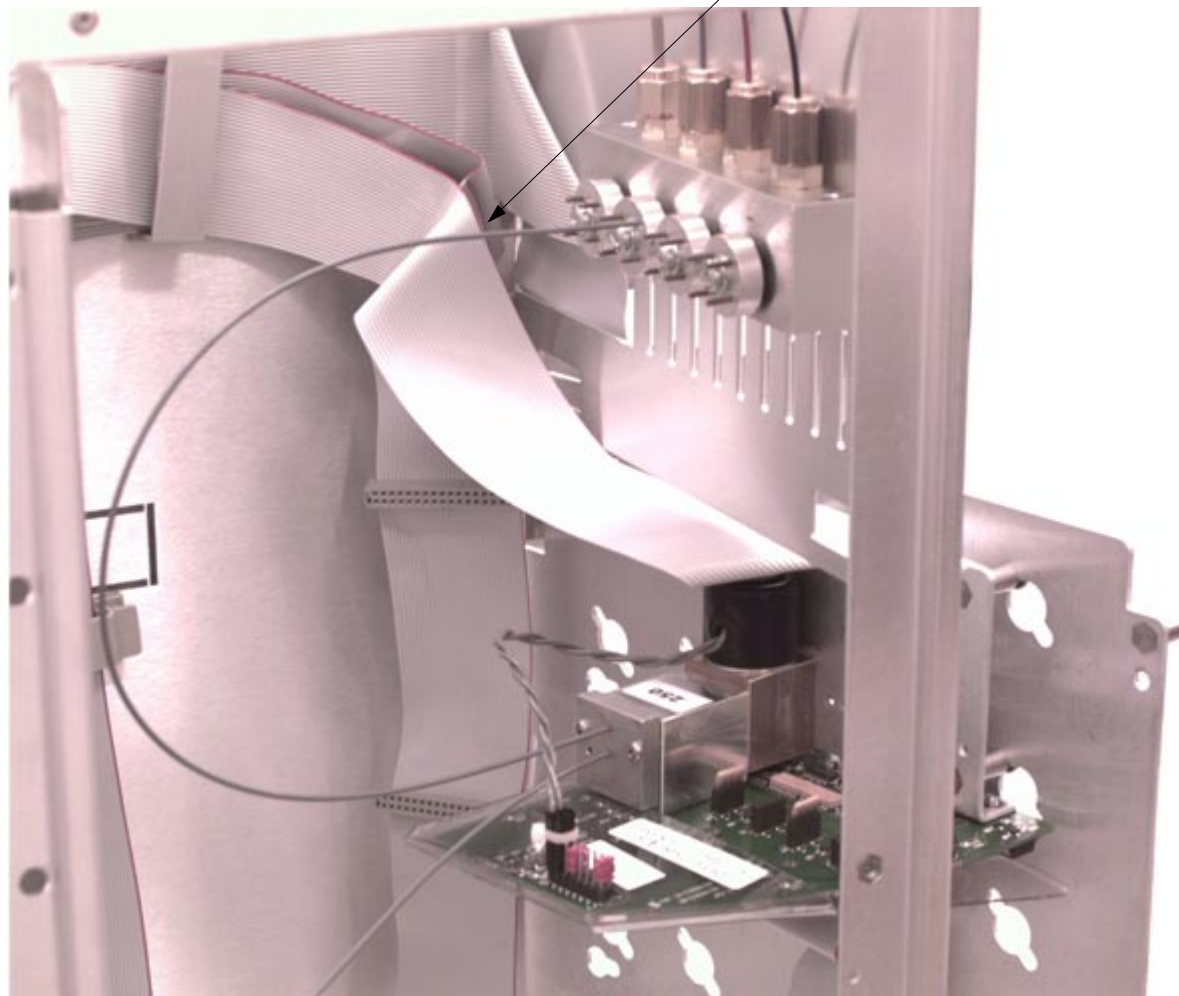

NOTES:

1. DRESS PLUMBING IN PNEUMATICS COMPARTMENT USING CABLE TIE. INSPECT TO ENSURE PNEUMATICS PANEL SWINGS FREELY.

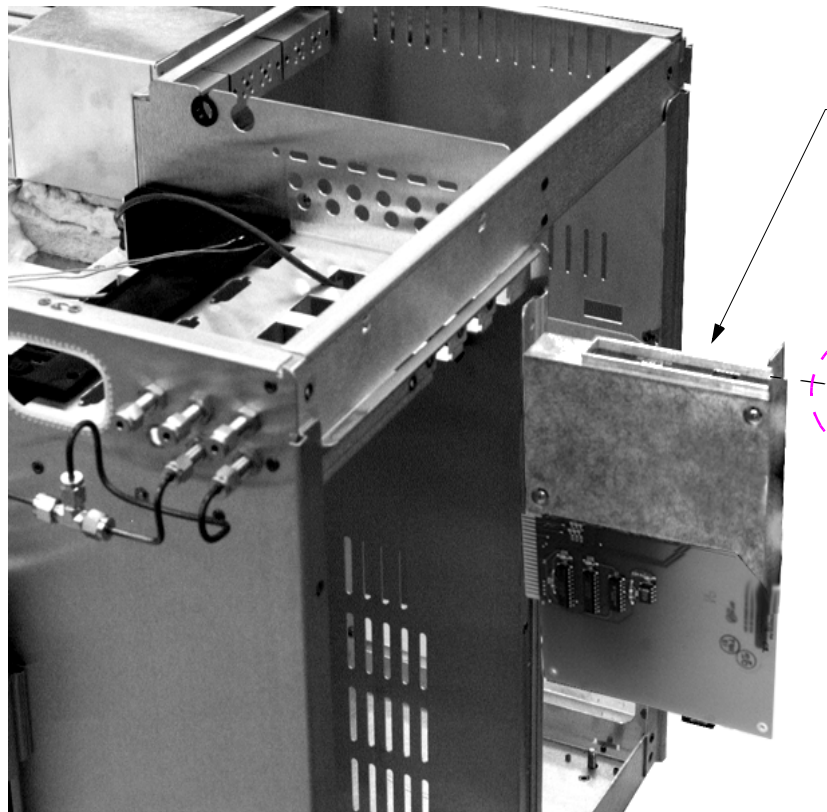

ECD PWA INSTALLATION


03-925021-01
FID PWA

03-925487-01
IGNITOR CABLE



03-917830-00
SIGNAL CABLE 18"

SIGNAL & IGNITOR CABLE CONNECTION

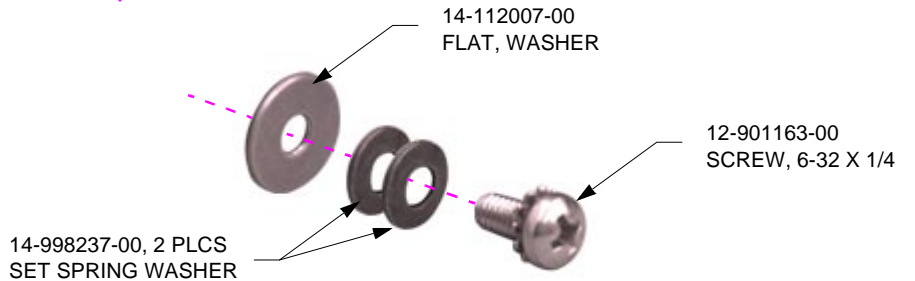

14-112007-00
FLAT, WASHER

12-901163-00
SCREW, 6-32 X 1/4

14-998237-00, 2 PLCS
SET SPRING WASHER

MOUNTING SCREWS DETAIL